Distributeur volume variable 10 mL

Réf: 451 005

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Read carefully before use!

Safety Instructions

This instrument is not designed for universal use. This instrument may be involved with hazardous materials, operations, and equipment.

This Manual does not purport to address all of the safety problems associated with its use. It is the responsibility of whomever uses this instrument to consult and establish appropriate safety and health practices and determine the applicability of regulatory limitations prior to use.

1. Every user must be acquainted with this Operating Manual before operation.
2. Observe the general instructions particularly for operating, cleaning and maintenance.
3. Follow the general instructions for hazard prevention and safety regulations, e.g., wear protective clothing, goggles and gloves.
4. Observe the reagent manufacturer's information and instructions.
5. Use the instrument only for dispensing liquids, with strict regard to the defined Operating Exclusions and Limitations (⇒ pages 6-7).

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If in doubt, contact the manufacturer or supplier.

6. When dispensing, the discharge tube must always point away from the user or any other person.
7. Attention! Due to the spring-operated mechanism, the dispensing cartridge will always be filled with liquid after initial priming.
8. Never carry the mounted instrument by the pump assembly or the valve block. Breakage or loosening of the plastic cylinder may lead to personal injury from chemicals.
9. Never depress the piston when the stopper cap is covering the discharge tube.
10. Never remove the discharge tube when the dispensing cartridge is filled with liquid.
11. Never use force on the instrument.
12. Use only original manufacturer's accessories and spare parts.
13. In case of trouble (e.g., piston difficult to move, sticking valves or leakage), immediately stop dispensing. Clean the instrument according to the cleaning instructions before any further use of the instrument (⇒ pages 14-16) or contact the manufacturer.

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Application

This instrument is designed for dispensing liquids, observing the following limits:

- 15 to 40 °C (59 °F to 104 °C) of instrument and reagent (other temperatures on request)
- vapor pressure up to 500 mbar
- viscosity:
25 ml instrument: 75 mm²/s
10 ml instrument: 150 mm²/s
- density up to 13.6 g/cm³

Valve block (N), filling tube (L), discharge tube (I) and valves (H, M) withstand steam sterilization at 121 °C.

The dispensing cartridge (B) and pump assembly (A) are not resistant to steam sterilization.

Operating range

- aqueous solutions of inorganic and organic substances
- alkaline solutions, e. g. ammonia, sodium hydroxide
- polar solvents, e. g. ethanol

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Please see our catalog for a listing of chemical compatibility with components of this instrument.

Please ask for further information.

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Operating Exclusions

Never use this instrument with liquids which attack PP, PE or EPDM.

Liquids that should not be dispensed include, but are not limited to:

- non-polar solvents like hydrocarbons and halogenated hydrocarbons
- concentrated or oxidizing acids.

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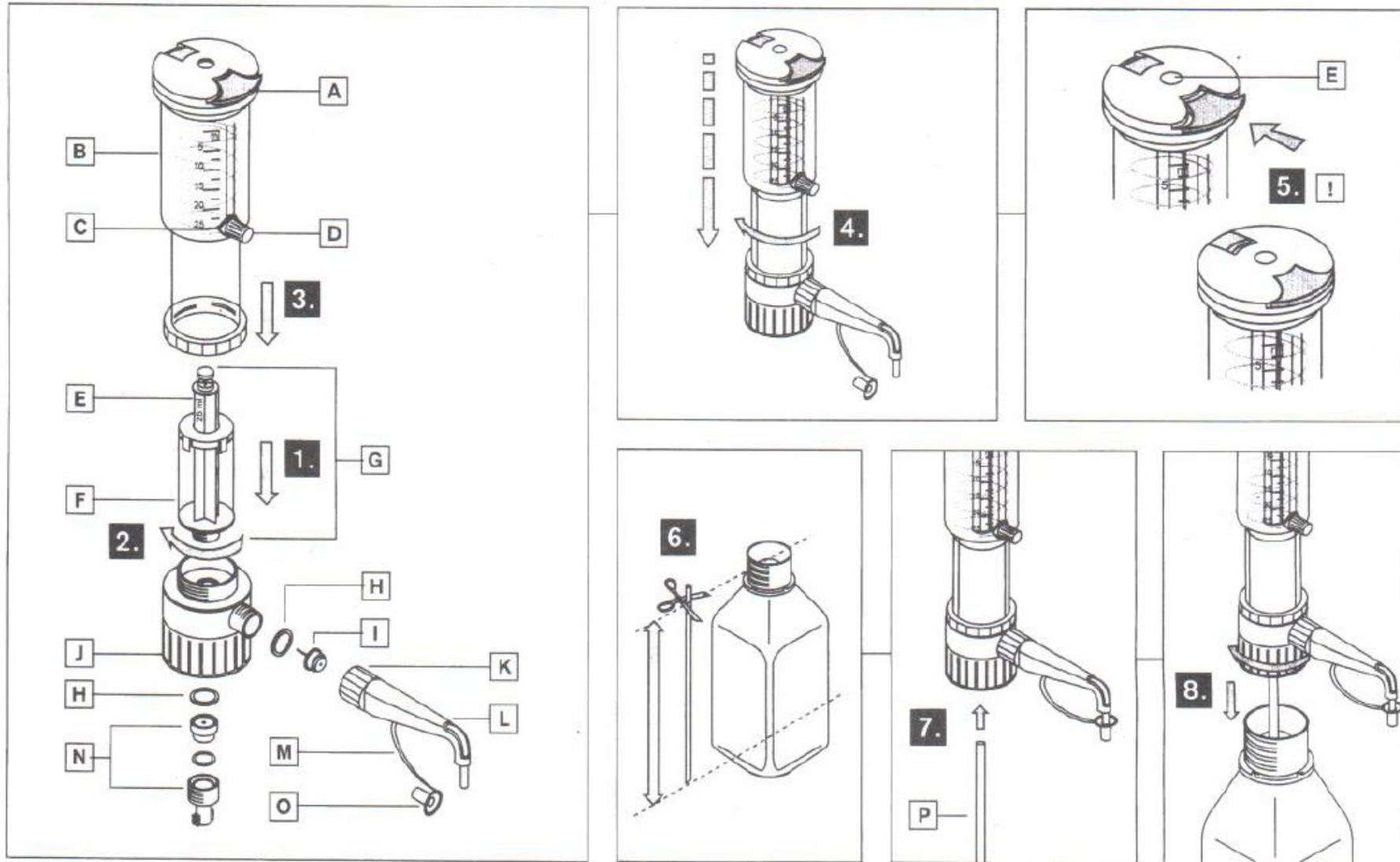
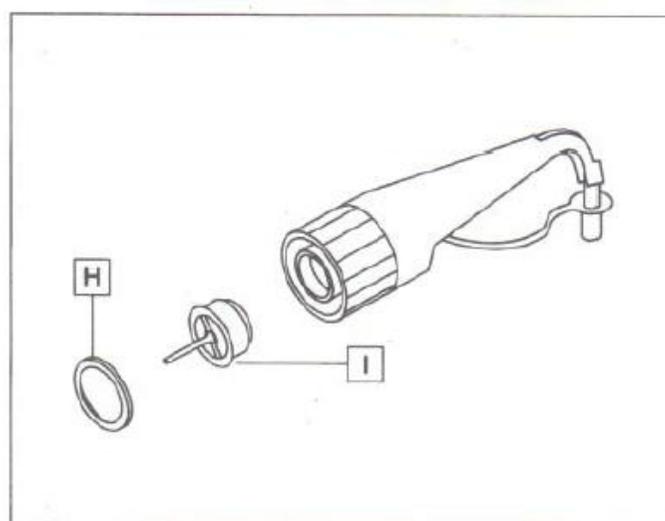
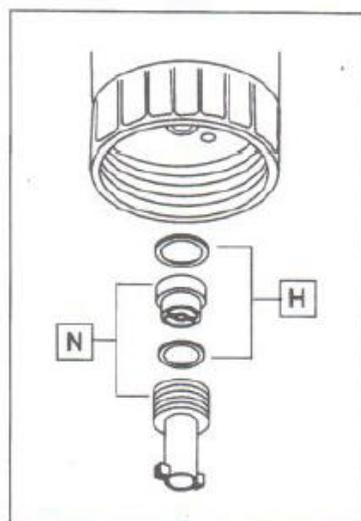
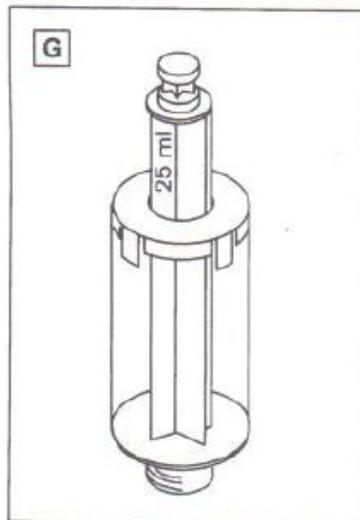
Components

- A Piston lock
- B Pump assembly, PC
- C Pointer
- D Volume adjustment knob
- E* Piston, PE**
- F* Dispensing cylinder, PP
- G* Dispensing cartridge, PP/PE
- H* Seals, EPDM
- I* Discharge valve, PP/EPDM
- J* Valve block, PP
- K Locking nut, PP
- L* Discharge tube, PP
- M* Stopper, PP
- N* Filling valve, PP/EPDM
- O* Stopper cap, PP
- P* Filling tube, PP

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Preparing for use

Warning!

Wear protective clothing, goggles and gloves. Follow all Safety Instructions and observe the Operating Exclusions (⇒ pages 4-7).

Note:

Follow all instructions on p. 8/9 for proper assembly.

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Priming the instrument and Volume setting

Attention!

Use a bottle stand when using smaller or other potentially unstable bottles to help protect against tipping. (Accessories and Spare Parts, ⇒ p. 23). Observe all Safety Instructions (⇒ p. 4-5).

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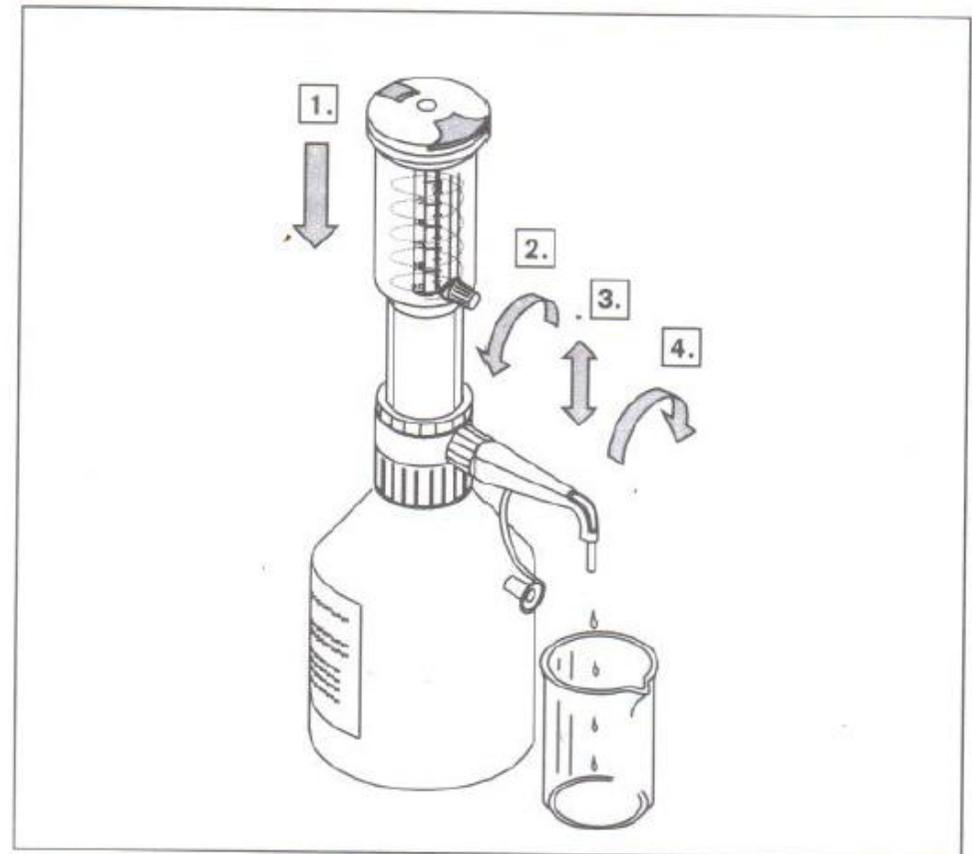
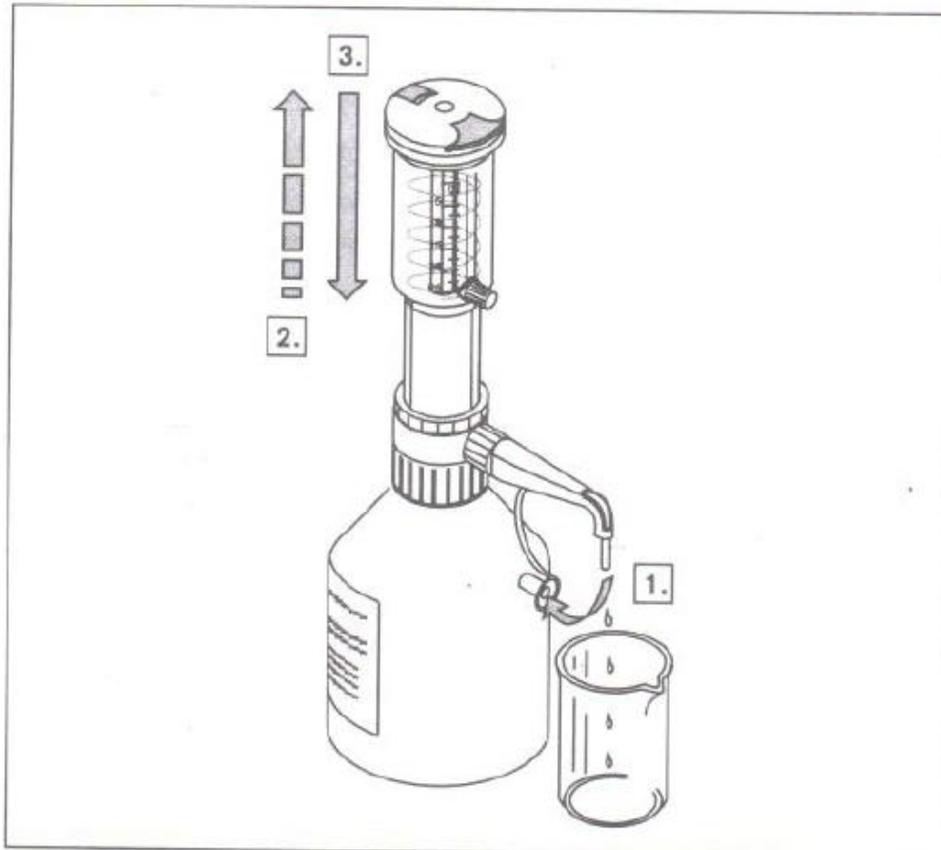
Dispensing

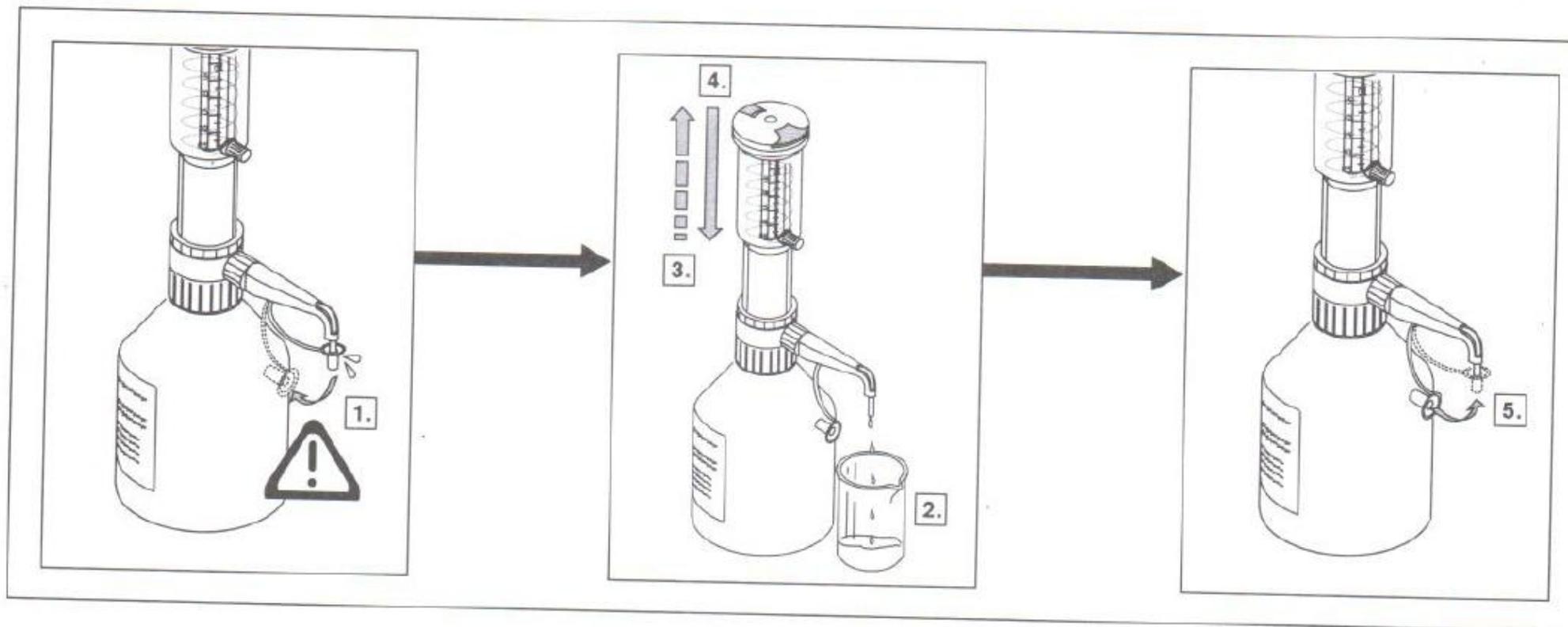
Attention!
Follow all Safety Instructions and
observe the Operating Exclusions
(⇒ pages 4-7).

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Replacing the dispensing unit

Attention!

The dispensing cartridge is a wearing part. Replace the cartridge when the piston is difficult to move or is not airtight. To prevent chemical injury, the dispensing cartridge has to be cleaned (⇒ pages 14-16) prior to replacement.

Note:

Follow the instructions on pages 8/9 for replacing the dispensing cartridge.

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Cleaning

The instrument must be cleaned in the following situations to assure correct operation:

- immediately if the piston becomes difficult to move
- before changing the reagent
- prior to long term storage
- before replacing the valves
- before replacing the dispensing unit

Depending upon the frequency of use, the dispensing cartridge should be replaced regularly.

Warning!

Follow all Safety Instructions (⇒ page 4-5).

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Warning!

Carry the mounted instrument only as shown in the figure.

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Note:

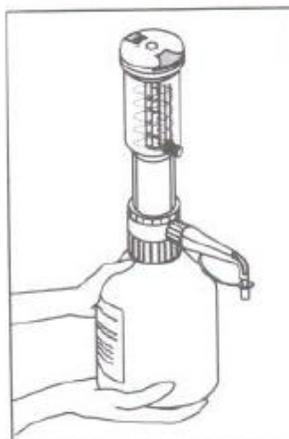
Clean only inside a suitable collecting basin.

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
- A** mild laboratory detergent, e.g., Biorex
- B** suitable solvent, e.g., distilled water

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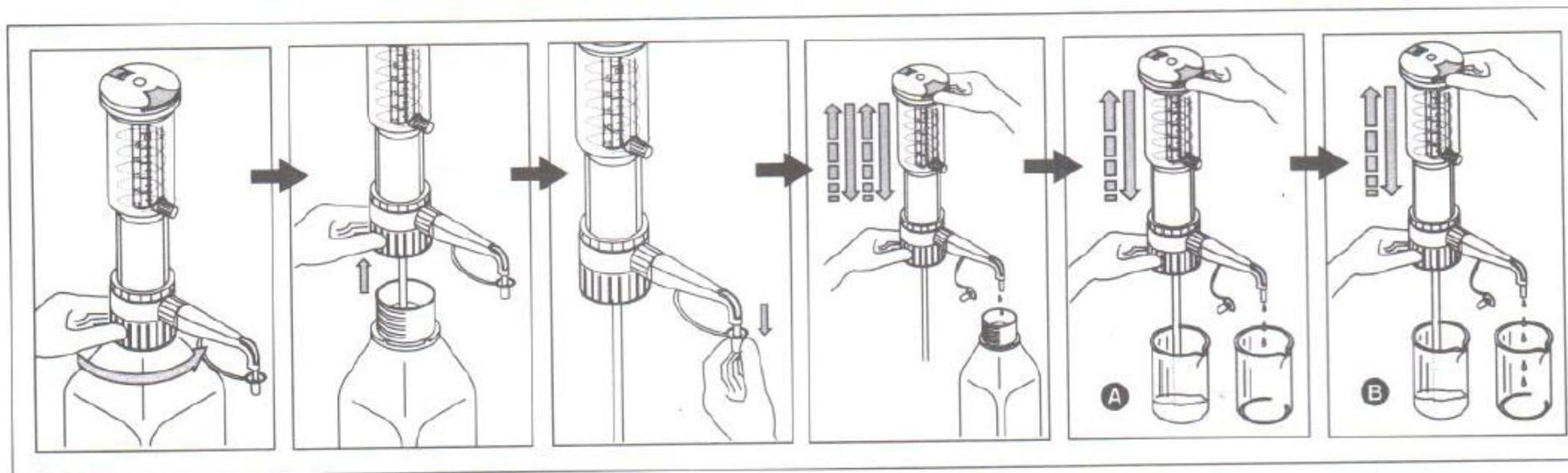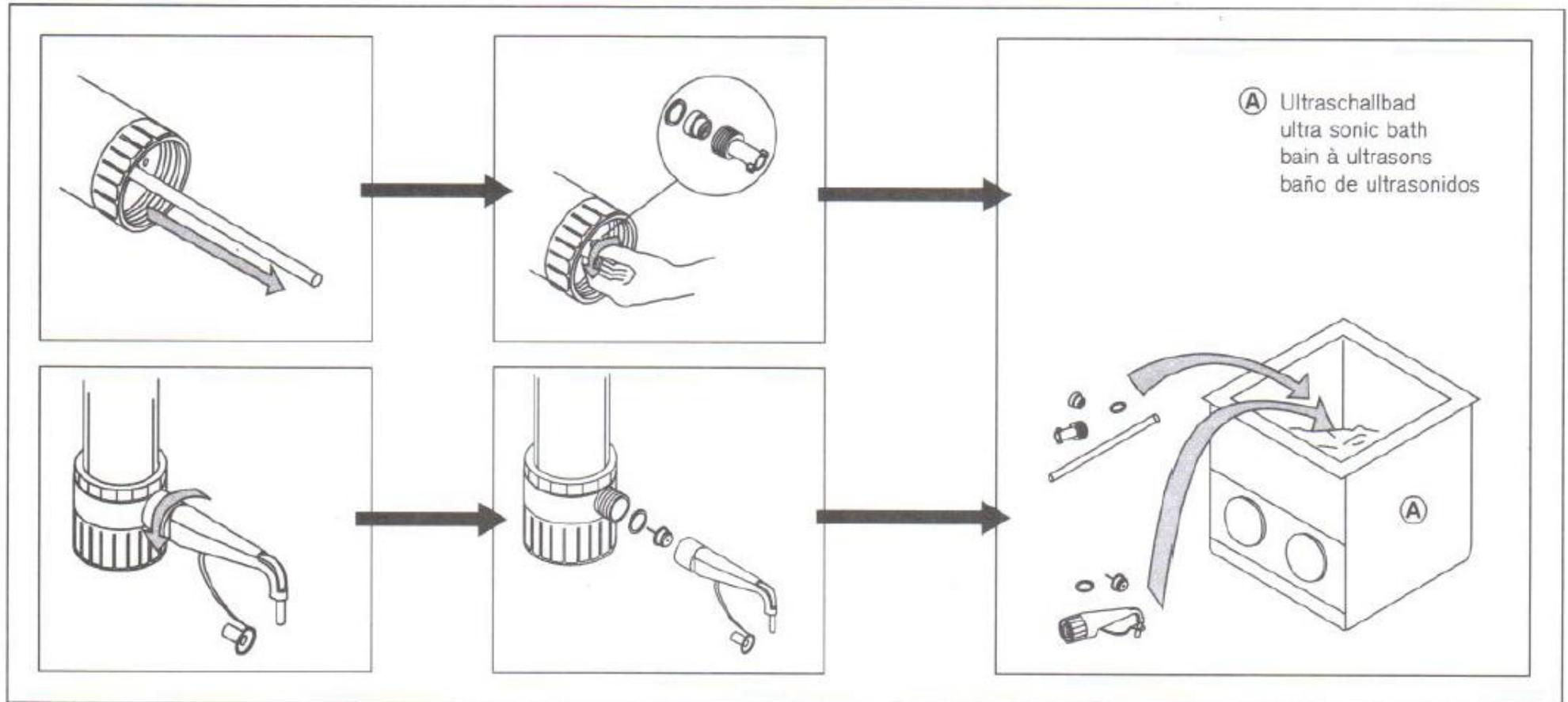
* The values of accuracy and coefficient of variation are final test values referring to the delivered volume, instrument and distilled water at equilibrium with ambient temperature (20 °C/ 60 °F) and smooth operation.
A = Accuracy
CV = Coefficient of variation

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